

Drinking Water Sample Set 1

IOCs - Inorganic Compounds
Metals Cyanide Fluoride

VOCs 524.2
Volatile Organic Compounds
No air in vials once they are filled.
Do Not Rinse! (Add 5 drops from the blue capped dropper to each vial)

Add contents of red capped vial.

Add contents of green capped vial.

Trip Blank
(Do Not Open)

No air in vials once they are filled. Do Not Rinse!

SOCs - Synthetic Organic Compounds

504	547	548.1	549.2	515.4	525.2	531.1
EDP, DBCP Pesticides	Herbicides	Semi-Volatile Compounds	Pesticides	Herbicides	Semi-Volatile Compounds	Carbomate Pesticides

Note: All of the above SOC bottles are pre-preserved, DO NOT RINSE!



No air in vials once they are filled. Do Not Rinse!

Add contents of blue capped vial to each bottle.

V. 12/14